## IDEXX SNAP\* Beta-Lactam ST Test

### Penicillin G detection level at or below 4 ppb

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**The SNAP\* Beta-Lactam ST Test is validated to detect residues in cow, goat, sheep and reconstituted whole fat powdered milk. Not for use in NCIMS milk-regulatory programs. Contains plastic, biological, and other components; in vitro use only. Not for sale in the U.S. Use test before expiration date.**

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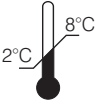
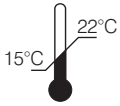
## Ensure test and sample are at correct temperature before use.

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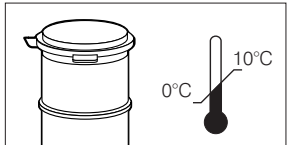
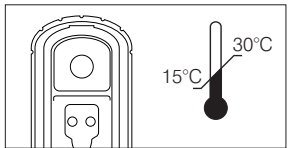
**Precautions:** Thoroughly mix milk before testing. Do not use abnormal-looking milk or expired tests; do not mix components from different kits. If the control spot fails to develop color, retest the sample.

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### 1 Transfer milk and swirl until dissolved, but no more than 15 seconds.

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

O  
OU  
OU  
LUB  
ODER  
OPPURE  
或  
または
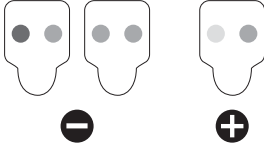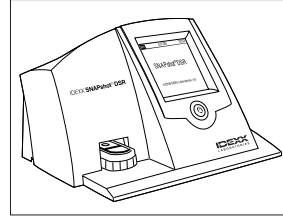
## IDEXX SNAPshot\* DSR Reader

### It is recommended that positive samples be further analyzed.

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### \*Stored at 15–22°C, test expires in 30 days or as labeled, whichever is earlier.

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### Technical assistance • [idexx.com/dairy](http://idexx.com/dairy)

US/Canada/Latinoamerica	+1 800 321 0207	日本	+021 6127 9528
EU	+00800 329 43399	Within AU	+1300 44 33 99
中国	+010 84423133	Within NZ	+0800 102 084

One IDEXX Drive • Westbrook, Maine 04092 USA

\*Samples of this test kit model were independently evaluated by the AOAC Research Institute and were found to perform to the producer's specifications as stated in the test kit's descriptive insert. The producer certifies that this kit conforms in all respects to the specifications originally evaluated by the AOAC Research Institute as detailed in the PERFORMANCE TESTED certificate number 061202.

Independently validated by  
**ILVO**

PERFORMANCE TESTED +  
**AOAC**  
RESEARCH INSTITUTE  
LICENSE NUMBER 061202

\*SNAP, SNAPshot and Simplicity from IDEXX are trademarks or registered trademarks of IDEXX Laboratories, Inc. or its affiliates in the United States and/or other countries.

Patent information: [idexx.com/patents](http://idexx.com/patents).

© 2013 IDEXX Laboratories, Inc. All rights reserved. • 06-19278-02

**IDEXX**